



587 East Middle Turnpike, P.O. Box 370, Manchester, CT 06040  
Telephone: 860.645.1102 Fax: 860.645.0823

## **Procedure for PCB Wipe Sampling**

*This technique can be used on hard, smooth, non-porous surfaces*

A wipe sample is taken by applying hexane to one gauze pad. The moistened gauze pad is then held with a pair of stainless steel forceps or rubber gloves and rubbed thoroughly over a 100 cm<sup>2</sup> area of the sample surface (delineated by a template) to obtain the sample. Place pad in the sample jar provided, which should be tightly capped and labeled.

If you have any questions, give us a call!

**860-645-1102**